


MEASUREMENT AND CONTROL

DIGITAL TEMPERATURE INDICATOR

These are generally used as headers of shell and tube heat exchangers and columns.

This instrument is mainly used to monitor the temperature of liquid in a glass vessel in a typical Glass Distillation Unit.

The instrument consists of a Temperature indicator and a Resistance Temperature Detectors (RTD). The instrument works on 230V, 50Hz power supply. This displays the temperature in degree Centigrades in three and half digits of 12.5mm character height.

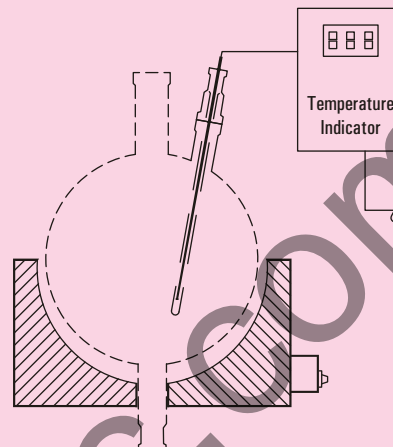


Cat.Ref.	Vessel size	RTD Length
DTI20	20	400
DTI50	50	500
DTI100	100	600
DTI200	200	700

TWO POINT DIGITAL TEMPERATURE INDICATOR

This instrument is mainly used to monitor the temperature of liquid in a glass vessel and temperature of vapours at reflux divider in atypical Glass Distillation Unit.

The instrument consists of a Temperature indicator and two Resistance Temperature Detectors (RTDs). The instrument works on 230V, 50Hz power supply. This displays the the temperature in degree centigrades in three and half digits of 12.5mm character height. A switch is provided to see the two temperatures alternatively.

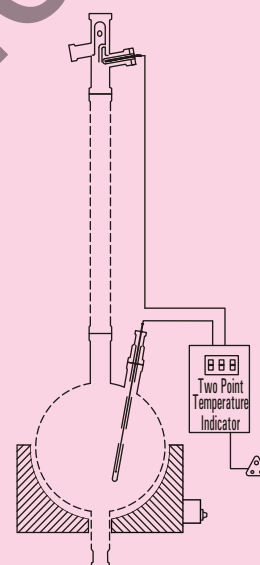


Cat.Ref. Length	Vessel size	RTD Length for vessel	RTD for reflux divider
DTT20	20	400	200
DTT50	50	500	225
DTT100	100	600	250
DTT200	200	700	300

ELECTRO-MAGNETS

Electro - magnets are used to operateMagnetically operated Reflux dividers. When 'On' the magnet attracts the swinging funnel of the reflux divider so that distillate can be taken off. Electro-magnets are to be mounted outside the glass column, just near to the reflux divider,with the help of adjustable fittings. These are designed to use with Timers to maintain correctratio between 'Off' and 'On' timings of its activation.

Electro-magnets work on 220V DC power supply,for which a output socket is provided in the Timers.

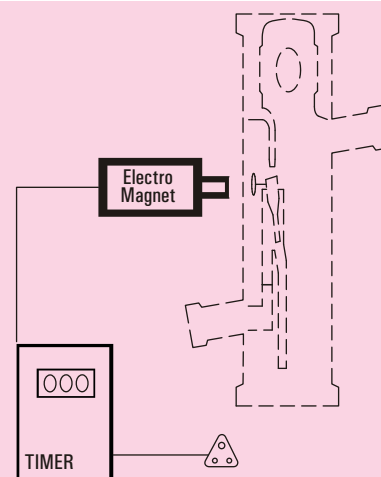

Cat. Ref.	Type
RPM*	Non-flameproof
RPF	Flameproof

TIMERS


Timers are designed to use with Electro-magnets to provide a correct ratio of refluxand distillate when operating a Magnetically operated reflux divider.

Timers work on a power supply of 230V, 50Hz.

Cat. Ref.	Type
RPM*	Non-flameproof
RPF	Flameproof

MEASUREMENT AND CONTROL

CONTINUOUS TEMPERATURE CONTROLLER

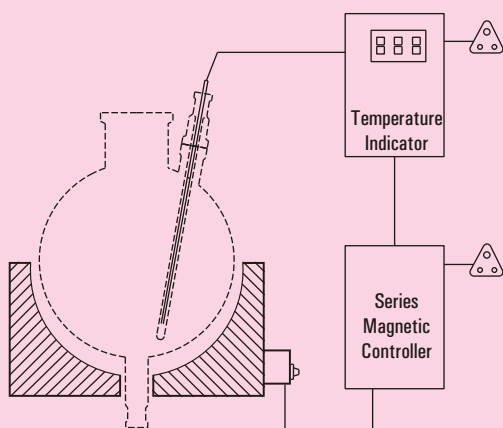


This instrument displays and controls the temperature continuously by switching the power supply on and off in an electrical heating equipment as per the initial settings of heating temperature, band width and reset temperature.

The instrument consists of a Temperature controller, a series magnetic controller & a Resistance Temperature Detector (RTD). RTD is put into the Thermometer Pocket of the glass vessel and desired settings are done. As the temperature in the vessel reaches to the set heating temperature, the temperature controller cuts the power in heating equipment off. Power starts on again as the temperature goes down as per the settings of band width and reset temperature.

This instrument works on power supply of 230V 50Hz and can be used with heating mantles and heating baths of all the size. It displays the temperature in degree centigrades in three and half digits of 12.5mm character height.

Cat.Ref.	Vessel size	L
CTC20	20	400
CTC50	50	500
CTC100	100	600
CTC200	200	700

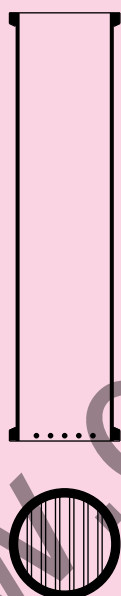

COLUMN SECTION WITH INBUILT PACKING SUPPORT

Goel introduce single piece column section with inbuilt packing support.

Advantages of this column against conventional column section:

- Ease in installation being a single piece instead of two pieces.
- There is no need to maintain stock of CS and packing support.
- Increases effective packed height which results in to increase in efficiency.
- Zero maintenance against column flooding.